



Spectradyne nCS1

Installing & Updating the Spectradyne Tools Software Package

© Spectradyne LLC

Vers. 2.0.1, 10/25/2016

Instructions for Installing/Upgrading the Spectradyne Tools Software Package

1. Uninstall prior versions

- To check the version and release date of the currently installed software, run:
Start > Spectradyne Tools > About
- Uninstall any old versions of Spectradyne Tools using:
Windows Settings > System > Add/remove programs

2. Download the software

- Download the Spectradyne Tools package from the following link:

<http://www.nanoparticleanalyzer.com/support.php>

3. Install Spectradyne Tools

- Run the setup.exe file to begin the install process
- Follow the steps in the installer

4. Update the calibration file on the installation computer

- Connect to the internet and run Spectradyne Tools > updateCal
- This App will download the latest available file called cartridgeCal_latest.csv to the location:
C:\Spectradyne\reffiles
- ** Note: If the computer is not connected to the internet:
 - o From a computer with internet access, download the file CartridgeCal_latest.csv from the Spectradyne website: www.nanoparticleanalyzer.com/support.php
 - o Copy the downloaded file to the following location on the installation computer:
C:\Spectradyne\reffiles
- This file is required to process data in the Spectradyne Tools Viewer App

5. Run the Software

- Use the desktop icons or start menu to start the Viewer or nCS1 software
- ** Note: The nCS1 software requires connection to an instrument before use

More Information and Technical Support

More detailed information on the operation of the nCS1 and its software can be obtained by sending email to support@SpectradyneLLC.com.

At Spectradyne we are eager to help with your particle analysis, so please contact us for help with specific measurement challenges and for help with the analysis of specific particle populations.



Speed. Precision. Range.

info@SpectradyneLLC.com

www.SpectradyneLLC.com